

Marine Broth (MB) medium

Into 100 mL distilled water dissolve 3.74 g powder medium of Difco™ Marine broth 2216 (Becton, Dickinson and Company) and 25.2 mg powder medium of Daigo IMK (Nihon Pharmaceutical Co., Ltd.). In the NIES-Collection, MB medium is used after autoclaving (121°C, 20 min).

Reference

Ishii, Y., Maruyama, S., Fujimura-Kamada, K., Kutsuna, N., Takahashi, S., Kawata, M., Minagawa, J. 2018 Isolation of uracil auxotroph mutants of coral symbiont alga for symbiosis studies. *Sci Rep*, **8**, 3237.