

Media for protozoa

Hemi

Into 98.9 mL distilled water 3.6 g powder medium of Daigo's Artificial Seawater SP (Nihon Pharmaceutical Co., Ltd.) and 0.1 mL LB broth medium (BD). After autoclaving, add aseptically 10 mL horse serum (Sigma-Aldrich Co. LLC).

Reference

Tashyreva, D., Prokopchuk, G., Yabuki, A., Kaur, B., Faktorová, D., Votýpka, J., Kusaka, C., Fujikura, K., Shiratori, T., Ishida, K., Horák, A., Lukeš, J. 2018 Phylogeny and morphology of new diplomonads from Japan. *Protist*, **169**, 158-179.