

## Media for protozoa

### **YE100ppm**

Into 100 mL distilled water 9 mg powder medium of Daigo's Artificial Seawater SP ( Nihon Pharmaceutical Co., Ltd. ) and 10 mg yeast extract.

## Media for protozoa

### P IV metals

Na <sub>2</sub> EDTA · 2H <sub>2</sub> O	100 mg
FeCl <sub>3</sub> · 6H <sub>2</sub> O	19.6 mg
MnCl <sub>2</sub> · 4H <sub>2</sub> O	3.6 mg
ZnCl <sub>2</sub> <sup>1)</sup>	1.04 mg
CoCl <sub>2</sub> · 6H <sub>2</sub> O	0.4 mg
Na <sub>2</sub> MoO <sub>4</sub> · 2H <sub>2</sub> O	0.25 mg
Distilled water	100 mL

1) In the NIES-Collection, 1.04 mg ZnCl<sub>2</sub> is replaced by 2.2mg ZnSO<sub>4</sub> · 7H<sub>2</sub>O.

### Reference

Provasoli, L., Pintner, I. J. 1959 Artificial media for fresh-water algae: problems and suggestions. In *The Ecology of Algae. Spec. Pub. No. 2.*, Eds. by Tryon, C. A., Jr. & Hartmann, R. T., Pymatuning Laboratory of Field Biology, University of Pittsburgh, Pittsburgh, p. 84-96.